

# INSTECH


## Covance Infusion Harness™ for Mice Model CIH62

Instech Laboratories, Inc.  
5209 Militia Hill Road, Plymouth Meeting, PA 19462-1216  
800-443-4227 • 610-941-0132 • FAX 610-941-0134  
EMAIL support@instechlabs.com • WEB www.instechlabs.com

### To assemble...

- Insert spring by twisting clockwise and pushing into top of saddle.\*
- Attach belly bands. To feed second end of belly band through center hole, wet tubing, push tubing end onto metal tool, and push through center hole as shown. The snug fit will hold the bands in place.\*
- Feed cannula up through tether, attach to swivel.

\* If preassembled, skip these first 2 steps.

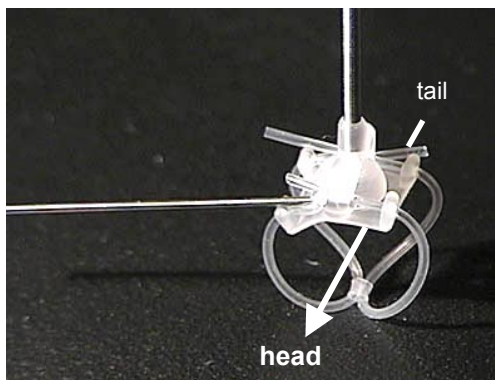

### To attach to the animal...

- Place loosely assembled harness around the animal. The front legs should fit between the belly bands and the spine should be oriented as shown above.
- Pull tube ends to tighten. Be sure harness is not too tight that it irritates the mouse, nor too loose that it moves around.
- Trim excess tubing, but leave some room for the animal to grow. Check tension regularly and adjust as needed.