


Model 400 Pump Mounting System

P400/MS Instructions for Use


INCLUDED COMPONENTS

- C-clamp with V-lock base station
- Aluminum pump mounting plate with rubber feet
- V-lock wedge
- (2) 1/4-20 pump mounting screws
- (3) 4-40 3/8in screws for V-lock wedge
- Screw driver
- 5/32 allen key

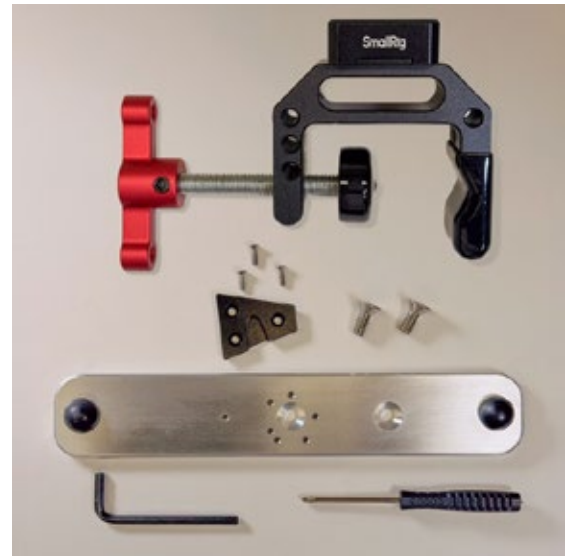

INSTALL MOUNT ON PUMP

1. Attach mounting plate to base of pump using large screws.

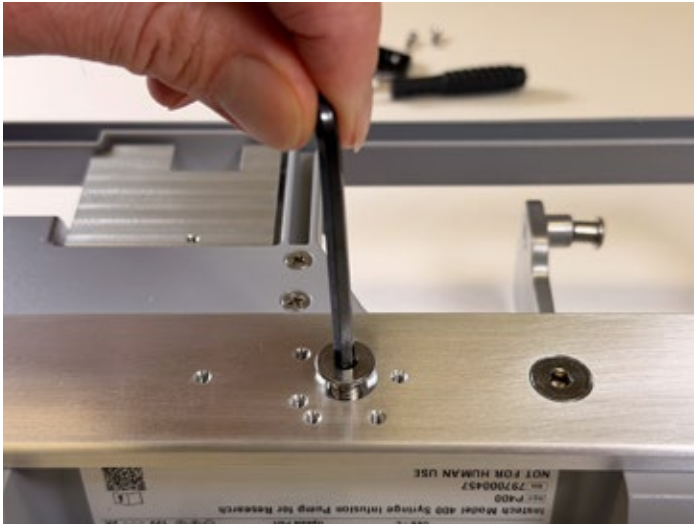


2. Attach V-lock wedge to mounting plate using small screws as shown to mount vertically on pole or horizontally on shelf. Use other mounting holes for other orientations.



INSTALL C-CLAMP

Pole Mount

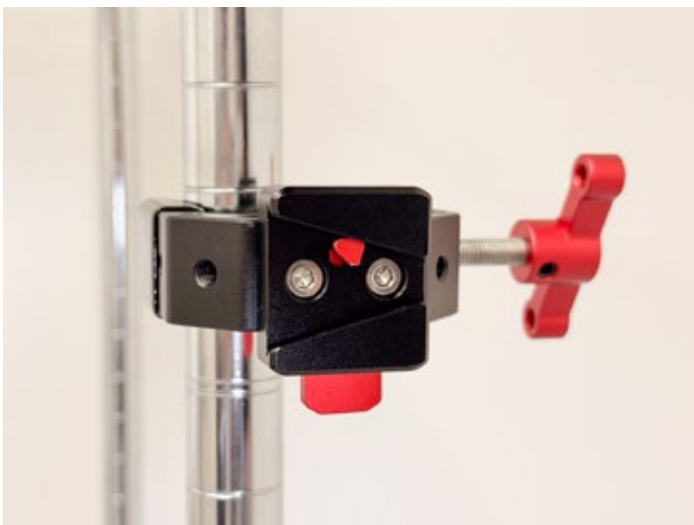

Shelf Mount



ENGAGE PUMP WITH C-CLAMP

Pole mount with vertical pump

(Note: never orient pump with syringe facing up.)



Shelf mount with horizontal pump



REMOVE PUMP FROM C-CLAMP

1. Locate red release button



2. While holding pump with other hand, press release button



INSTECH

Instech Laboratories, Inc. | Plymouth Meeting, PA USA | 800.443.4227 | 610.941.0132
www.instechlabs.com

Part number: 7700-05-0216